

Calculate ratiometric FRET-Images

Summary

This tutorial shows step-by-step, how the "FRET Image" script of SymPhoTime 64 can be used to calculate pixelby-pixel the average FRET efficiency in an image containing several cells transfected with Cerulean/YFP constructs. To calculate the average FRET efficiency, a ratiometric approach based on the intensities in the donor and the acceptor channel is used.

Background Information

FRET is a method to determine molecular interactions or conformational changes based on a suited donor/acceptor pair.

Intensity FRET as used here can only be used to calculate average FRET efficiencies. In case of a mixed population, these populations can only be separated using a donor dye with a single exponential decay kinetic and analyzed with lifetime data. This is not the scope of this tutorial.

The script usage is demonstrated on an image of several cells which have been transfected with CENP-B-Cerulean and YFP-CENP-A. Cerulean is a CFP variant an frequently used as a FRET donor. The CENP-complex consists on several proteins and is involved in mitosis. For further details on the sample, see S. Orthaus et al., Chembiochem. 2008, 9, 77-92, and the PicoQuant application note

http://www.picoquant.com/images/uploads/page/ files/7267/appnote_flim_fret.pdf.

The example image shows four cells, two of them were only transfected with CENP-B-Cerulean, the other two contain both constructs. The same image is also used for the tutorial of the lifetime FRET image script. Check there for comparison.

In ratiometric FRET, a lot of system parameters have to be determined from several images. This problem is too complex to be accessed with a single tutorial. In this tutorial, some arbitrary correction parameters are used for demonstration purposes only.

Several approaches have been published as procedures to calculate the needed correction parameters, e.g. ¹; ². Evaluating and reviewing all papers in this area would overpass the scope of this tutorial.

Step-by-Step Tutorial

- Start SymPhoTime 64 software.
- Open the "Samples" workspace via "File\open Workspace" from the main menu.

Note: The "Samples" workspace is delivered with the SymPhoTime 64 and on the DVD-ROM and contains example data to show the function of the SymPhoTime 64 data analysis. If you haven't installed it on your computer, copy it from the DVD onto a local drive before going through this tutorial.

Response: The files of the sample workspace are displayed in the workspace panel on the left side of the main window.



• Highlight the file CENP-labelled _cells_for_FRET.ptu by a single mouse click.



• Select the "Analysis" tab and in there, open the drop down menu "Imaging".

	Analysis	
	Imaging	
B-all Cy5_diff_RF+FLCS-pattern.ptu		Multi France Fi 114
⊕ Atto655_diff_FLCS-pattern.ptu	FLIM	Multi Frame FLIM
Atto655+Cy5_diff_FCS+FLCS.ptu	A FLIM analysis. Select an	A FLM analysis of individula
III Atto855_diff_2FFCS.ptu	image TITR file and press	frames. Select an image
Atto488_diff_cw_total_correlatio	Start.	TTTR file and press Start.
Atto488_diff_pulsed_antibunchin	C Halo L Stat	Hala I Start
Cy5_immo_FLM+Poi-imaging.ptu	Help Start	neip start
E-dd Cy5_immo_Lifetime_Trace.ptu	Grouped FLIM	Anisotropy Image
H-ALA Altoess_mmo_On-Off-Analysis.p	A Grouped FLIM analysis.	An intensity based
DaisyPollen_cells_FLM.ptu	Select any collection of	anisotropy image. Select a
GPP_RPP_cells_FLM-FRET.ptu	image TITR files and press	two channel image TTTR file
TS-Bead immo xy-scan Dual Fo	Tield	and press Start.
TS-Bead immo xz-scan.ptu	nep Start	neip Start
BA488+BA547 unlinked mix.ptu	FRET Image	Lifetime FRET Image
BA488+547_crosslinked.ptu	An intensity based FRET	A lifetime based FRET image.
Atto488_485nm_pulsed.ptu	image. Select a two channel	Select an image TTTR file
GUVs.ptu	Starf	and press Start.
NV-Center_for_Antibunching_1.p	State	
INV-Center_for_Antibunching_2.p	Help Start	Help Start
- Att NV-Center_for_Antibunching_se	Pattern Matching	Position in Image
CENP-labelled_cells_for_FRET.ptl	A TCSPC pattern based	Mark the positions of
-alla CENP-labelled_cells_for_FRET_IR	decomposition. Select an	measured files in the
L CENP-labelled_cells_tor_PRET_IR	Starf	overview mage.
FRET_GFP and mRFP.ptu	Male V Made	Hala I Start
	Help Start	Help Start
	Time Trace	
	TCSPC	
	 FCS 	
	Grouped Analysis	
	 Alignment 	
	User Defined Scripts	

Note: the drop down menu can be opened and closed by clicking on the grey button on the left side of the header of the drop down menu:

• Start the "FRET Image" script by clicking on "Start".

FRET Image		
An intensity based FRET		
image. Select a two channel		
image TTTR file and press		
'Start'.		
Help Start		

Response: The FRET image script is applied to the file CENP-labelled _cells_for_FRET.ptu. Thereby, a new Window opens:

😥 lonPluTime M. (FMT image - CIM	(algebra, etc.) WE give				_ICX
20 File Edit View Settings Scripts A	nalysis Window Help				_ # ×
20 Tel Line Land, Salanga Graph & Graph S. 20 Tel Line Land Colling Control of Salary Sciences 21 Advances of the Salary Sciences of Salary Sciences 21 Advances of Salary Sciences of Salary Science	Available Cas 192 Cas 192	<u>20 µm</u>		80 µm	
Total Strenger: T			L/MT W1: 6.0 % % W2: 19.0 % % Stagl Rei 61 % Read/abb million 6 Address million 6 Address million 6 Stagl Rei 6 Stagl Rei	200 700 700 700 700 700 700 700 700 700	No. Constant of the second sec

Note: The window contains five different regions:

- 1. Upper left: imaging calculation analysis options.
- 2. **Upper center:** Fast FLIM image, displayed in false color scale. As the donor and acceptor channels have not been assigned yet, this image area is still empty.
- 3. **Upper right:** FRET image. As it has not been calculated yet, this area is still empty.
- 4. Lower center: Here the FRET efficiency histogram over the selected pixels will be plotted when the calculation is performed.
- 5. Lower right: Here the FRET distance histogram in terms of R0 will be plotted when the calculation is performed.
- Set the pixel "Binning" to 2 points.
- As data channels, check only Channel 1 and 2, as in these channels the relevant photons are recorded.



• Press "Calculate".

Response:

• Automatically, the two active detector channels are assigned to the donor and acceptor channel. By default, detector channel 2 is considered the donor channel and detector channel 1 is assigned as acceptor channel.

Note: If this assignment is not correct, it can easily be inverted using the controls in the "Donor/Acceptor Assignment" panel.

• A false color image is calculated, in which the intensity of detector 1 and 2 are plotted in green and red,

respectively.

• Invert the color assignment to display the acceptor channel in red and the donor channel in green. Therefore, click on the bar next to the image to open the display options and invert the channel selection.



↓



Response: The image is adapted, the acceptor is now plotted in red.

Note: Now four cells are visible, the cells on the left are mainly plotted in green. These cells have only been transfected with the donor construct, while the other two cells are transfected with both constructs. The fluorescent spots are restricted to the nuclei, only some vesicles in the cytoplasm can be seen in the cell on the lower right.

- Set the Min. Photons per FRET event to "700" Cnts (counts).
- Click on "more" to open the additional options for entering the correction parameters.

 FRET Calc. 	Antion		
Min. Photons p	er FRET Event.		700 Cnts 🚔
	Calo	FRET	Help
(more)			
FRET Calculati	on		
Min. Dhotopo per E	DET Event		700 Cnts
ant, Photona per r	The F L F DOWN.		
ann, Photonis per r	Calc.	FRET	Help
ant, Photone per r	Calc.	FRET	Help
	Calc	FRET	Help
(less) Gamma:	Calc.	FRET	Help
Gamma:	Calc.	FRET	0,000
(less) Gamma: [Cross Taik: [Donor Background	Calc.	FRET	0,000

• As parameters, enter "70" for "Gamma" and "7" for "Cross Talk".

Note: For more information about the equation used for image calculation, use the help button next to the "Calc. FRET" button. This opens a window with the equations and a description of the fitting parameters. The correction parameters here are arbitrarily set to result in an image where the FRET efficiency matches roughly the values obtained in the "Lifetime FRET Image" script (see the corresponding tutorial). They have not been correctly determined and are only used for demonstrating the function of the script.

• Click "Calculate FRET".

↓

FRET Calculation					
	700 Cnts				
Calc. FRET	Help				
00					
00 Direct Exc.:	0,000 💂				
	0,00 Cnts				
	0,00 Cnts				
	Calc. FRET				

Response:

- The FRET image is calculated, and the corresponding histograms are calculated. In the FRET image the FRET efficiency is only calculated for pixels with intensities higher than the threshold.
- The FRET efficiency color scale is automatically set from 0 to 100% FRET efficiency.
- The FRET efficiency histogram and the Distances histogram are shown.



• It can be seen that a few fluorescent dots on the lower right are over the set intensity threshold. As these signals are not in an area of interest, place the mouse cursor over the FRET image, activate the context menu with a right mouse click and select the "Rectangle ROI" tool.



• Now it is possible to remove the unwanted area by constantly pressing the "CRTL" button and the left mouse button, draw a rectangle over these areas and release mouse and keyboard. Do this repetitively until all unwanted signals are removed.



• Press "Calculate FRET" again to update the histograms.

Response: The FRET efficiency and distance histograms are updated. This can be seen clearly at the small peak around 20% FRET efficiency which disappears now.

🔅 landaline M., PMT image - CM	National Society (Million)				_ICX
R File Edit View Settings Scripts A	nalysis Window Help				_ a ×
20 Service 20 For 16th Control Level 1. (MI Transp. Control Level 1. Control Level 1. Conttore 1. Co	Access Function Form (See) (Mit Cale) Access Function Form Access Function Form	100 µm		80 um	
Table Transport Torinstance Ground Angel Torinstanc		0.30 0.30 0.30 0.30 0.40 0.40 0.40 0.40	1,107 10. 10 % Her 10.1% Bel Re: 41 6 400000 hotigen 6 400000 hotigen 6 1000 Her 1000 1000 Her 1000 Her 1000 Her 1000 Her 1000 Her 1000 Her 1000 H		Deserve.

• Store the result by clicking on "Save Result.

🔺 File			
	Save	Result	
Sav	e Defaults	Restore Defaults	

Response: A result file is generated and linked to the raw data file.



- The image analysis is now finished. There are several possibilities to continue:
 - To adapt the histogram plots, change the plot parameters on the right of the histogram to the desired values and click "Recalculate". To save these plots, click "Save Results" again.
 - To export certain analyses, place the mouse cursor over the area of interest, activate the context menu with a right mouse click and choose the export function of interest to export the data.
 - In case one wants to analyze many images with the same FRET pair, it is possible to save the starting values for the fits by clicking on "Save Defaults" and store the starting values as a User Profile by entering the main menu ("Settings\Save User Profile as"). The software will need to restart to adapt the changes.
- If you don't want to do anything more, just close the result window.

¹⁾ Roszik et al., Cytometry A 2009, 75A, 761-767 ²⁾ Tadross et al., J. Microsc. 2009, 233, 192-204

Copyright of this document belongs to PicoQuant GmbH. No parts of it may be reproduced, translated or transferred to third parties without written permission of PicoQuant GmbH. All information given here is reliable to our best knowledge. However, no responsibility is assumed for possible inaccuraciesor omissions. Specifi cations and external appearances are subject to change without notice.



PicoQuant GmbH Rudower Chaussee 29 (IGZ) 12489 Berlin Germany P +49-(0)30-1208820-89 F +49-(0)30-1208820-90 info@picoquant.com www.picoquant.com