

Intensity Time Trace Analysis

Summary

This tutorial shows step-by-step, how to draw intensity time traces from a single molecule. The script requires a point measurement of an immobilized fluorophore. When applied to images, it can be used to check for line and frame marker presence and assignment.

Step-by-Step Tutorial

Select a file and start the script

- Start SymPhoTime 64 software.
- Open the "Samples" workspace via "File\open Workspace" from the main menu.

Note: The "Samples" workspace is delivered with the SymPhoTime 64 and on the CD-ROM and contains example data to show the function of the SymPhoTime data analysis. If you haven't installed it on your computer, copy it from the DVD onto a local drive before going through this tutorial, e.g. to a folder C:\SPT_User.

Response: The files of the sample workspace are displayed in the workspace panel on the left side of the main window.

SymPhoTime 64	
File Edit View Settings	Scripts Analysis Window Help
	Analysis
Samples	S-pattern ntu
in him cys_uni_ind free in him cys_uni_ind free in him Atto655_diff_FLC	S-pattern.ptu
tto655_diff_2FF(CS.ptu
_Atto488_diff_cw لىلىت Atto488_diff_cw	antibunching.pt
Cy5_immo_FLIM+F فالسلس Cy5_immo_Lifetim	Pol-Imaging.ptu e_Trace.ptu Grouped Analysis
Atto655_immo_On	-Off-Analysis. FLIM.ptu
GFP_RFP_cells_FI	IM-FRET.ptu
TS-Bead_immo_xy	/-scan_Dual Fo
TS-Bead_immo_xz	z-scan.ptu

• Highlight the file Cy5_immo_Lifetime_Trace.ptu by a single mouse click.

Note: This file contains a fluorescence time trace of an immobilized Cy5 molecule, its fluorescence was split by polarization onto two detectors.



• Select the "Analysis" tab and in there, open the drop down menu "TimeTrace".



Note: The drop down menu can be opened and closed by clicking on the grey button on the left side of the header of the drop down menu:

• Start the "Intensity TimeTrace" script by clicking on "Start".

由

¢

di la

Intensity Time Trace	
An intensity time trace. Select a TTTR file and press 'Start'.	
Help Start	

Response: The "Intensity Time Trace" script is applied to the file Cy5_immo_Lifetime_Trace.ptu. Thereby, a new Window opens:



Note: The window contains two different regions:

- 1. Left: Plotting options.
- 2. Center/right: Intensity time trace. As no time traces have been calculated yet, this area is still empty.
- Select a "Binning" of 10 ms.
- Click "Calculate".

 Trace Settings 			
Binning:			10,00 ms 🚔
Set Time Gate:			~1
, j	 		
Select Data Channels:			
1: 💌	2: 💌	3:	4:
	Calculate		

Response: As activated under "Select data channels", two intensity plots are calculated and plotted according to the two selected detector channels.



Note: The plot consists of several parts:

In the **small upper panel**, the complete time trace is plotted as the sum of all intensity channels. The **main intensity** time trace plots show the subsection of the trace marked in green in the upper trace for the two channels.

On the right, photon counting histograms (PCH) are plotted, indicating the frequency of different intensity values.

• To export data, e.g. the upper PCH histogram, place the mouse cursor in that region and open the context menu with a right mouse click. Select "Export ASCII: Selected cell".



Response: A window opens asking for a file name to store the data.

• Save the data. They can be opened e.g. with the "Notepad" program and imported into a suited fitting software.

D Export ASCII			1011	×
🕞 🕞 - 📧 Desktop		🔹 🏘 Deski	op durchsuchen	P
Organisieren 👻 Neu	er Ordner		≡ •	
 ★ Favoriten ■ Desktop Bownloads 32 Zuletzt besucht ⇒ Bibliotheken ⇒ Bider > Dokumente ↓ Musik ➡ Videos 	E Name	Große	Eementtyp	Änder
1 Computer	* K			
Dateigame: PCH	(CH1			•
Dateityp: ASC	I-Files (*.dat)			•
Ordner ausblenden		Spei	chern Abbre	chen

• Save the generated time traces by clicking "Save Result".

🔺 File		
	Save	e Result
	Save Defaults	Restore Defaults

Response: A result file TimeTrace.pqres is generated.

Samples
Ey5_diff_IRF+FLCS-pattern.ptu سلاس
Atto655_diff_FLCS-pattern.ptu السلاية
Atto655+Cy5_diff_FCS+FLCS.ptu فللس
Atto655_diff_2FFCS.ptu فللس
Atto488_diff_cw_total_correlation.ptu الملد
Atto488_diff_pulsed_antibunching.ptu الملك
Cy5_immo_FLIM+PoI-Imaging.ptu
Cy5_immo_Lifetime_Trace.ptu الملك
Cy5_immo_Lifetime_Trace.pco
Blinking_TimeTrace.pqres
Blinking_LifetimeTrace.pqres
TimeTrace.pqres
Atto655_immo_On-Off-Analysis.ptu المله
DaisyPollen_cells_FLIM.ptu
GFP_RFP_cells_FLIM-FRET.ptu
🖳 سلله Cy3+Cy5_diff_PIE-FRET.ptu
TS-Bead_immo_xy-scan_Dual Focus.
i → TS-Bead_immo_xz-scan.ptu

• The intensity time trace window can now be closed.

Checking Marker Signals in Images

Note: PicoQuant photon counting boards can register up to four external marker signals. These marker signals are usually generated by a scanning device. Usually, in a FLIM&FCS upgrade kit for Laser Scanning Microscopes (LSM), three markers are used: "line start", "line stop" and "frame marker".

These markers can be visualized using the "Intensity Time Trace" script to assign them correctly during the configuration of new setups.

• In the sample workspace, select the file GUVs.ptu and apply the time trace script to that file. When the intensity time trace window opens, click "Calculate" using a binning of 1 ms as set by default.



Note: When the intensity time trace of an image is displayed, the scanning results in intensity spikes.

• Open the pulldown menu "Marker" and select all markers.

Markers	
Marker 1 (51204)	
Marker 2 (51204)	<u> </u>
Marker 3 (101)	
Marker 4 (0)	

Response: The marker signals are drawn in the time trace. It can already be seen that the blue marker is more sparse. The marker pull down window also shows, now many markers are present.

Note: The image contains three different markers. Marker 1 and 2 are the most frequent markers, i.e. they serve as line start and line stop signals, whereas marker 3 is less frequent and serves as frame marker. In the image below, the frame markers can be clearly distinguished, while the line markers are too closely spaced for discrimination.



Zoom into the time trace image to distinguish between the individual line markers. This is done by
moving the mouse cursor into the region to be zoomed in, pressing the left mouse button and keeping it
pressed while moving to the cursor to the end of the region to be zoomed in and releasing the mouse
button. This has to be done several times until a sufficient magnification as in the image below is
achieved.



Note: It can be clearly seen that the distance between the black marker 1 and the next red marker 2 to the right is bigger than the distance between the red marker 2 and the following black marker 1 on the right. As this was a mono-directionally scanned image, the return from the end of the line to the line start position is done with highest possible speed. In contrast, the scanning of the actual line is performed according to the user settings, which are usually slower, but never faster than the backfly of the scanner. Therefore, in this measurement, marker 1 and 2 are the line start and line stop markers, respectively.

• You can zoom out again by activating the context menu with a right mouse click and select "Rescale".

• In this way, the marker assignment for this file has been checked and the window can be closed. As we are not interested in the data itself, there is no need to store this file.

Copyright of this document belongs to PicoQuant GmbH. No parts of it may be reproduced, translated or transferred to third parties without written permission of PicoQuant GmbH. All information given here is reliable to our best knowledge. However, no responsibility is assumed for possible inaccuraciesor omissions. Specifi cations and external appearances are subject to change without notice.



PicoQuant GmbH Rudower Chaussee 29 (IGZ) 12489 Berlin Germany P +49-(0)30-1208820-89 F +49-(0)30-1208820-90 info@picoquant.com www.picoquant.com