

~~TOC~~

## Phasor Analysis


Phasor Analysis is currently not included in [SymPhoTime64](#). However, you can use [Globals](#) developed by the [Laboratory for Fluorescence Dynamics](#) to analyse FLIM data via the Phasor approach.

## Install SimFCS

- Download [Globals for Images](#)
- Install SimFCS according to the documentation: <http://www.lfd.uci.edu/globals/>

## Exporting from SymPhoTime64

- Start FLIM analysis
- Calculate a fastFLIM image
- Select **Export Binary** <sup>1)</sup>

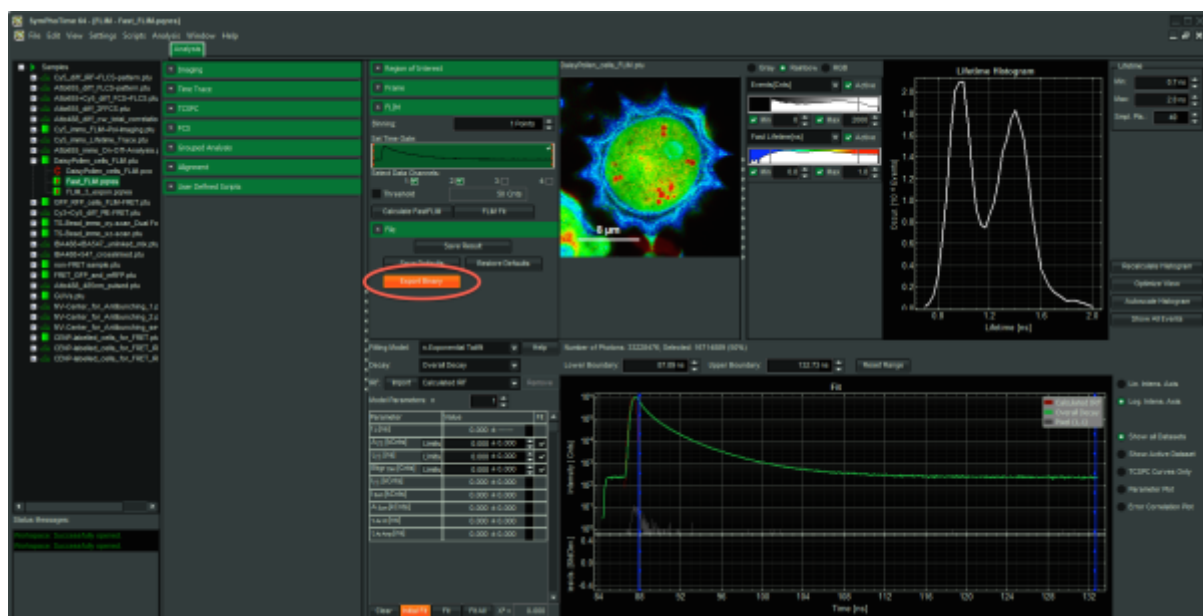

## Analysis in Globals/SimFCS

- follow the [Reading PicoQuant BIN files](#) tutorial of the Globals Software package to import the BIN files.
- follow the [Phasor Analysis Tutorial](#) of the Globals Software.
- [Deriving the laser frequency from PicoQuant BIN files](#)
- [Function: Read the PicoQuant file](#)

Users of SymPhoTime 32 with a PicoHarp 300 ,please refer to:

- [Using PicoQuant PT3 format for FLIM](#)
- [Reading PicoQuant PT3 files for FCS and lifetime FCS](#)

Additional Tutorials for the Globals software can be found here: <http://www.lfd.uci.edu/globals/tutorials/>

1)

## [Structure of the pre-histogrammed Image Data File](#)

Copyright of this document belongs to PicoQuant GmbH. No parts of it may be reproduced, translated or transferred to third parties without written permission of PicoQuant GmbH. All information given here is reliable to our best knowledge. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications and external appearances are subject to change without notice.



PicoQuant GmbH  
Rudower Chaussee 29 (IGZ)  
12489 Berlin  
Germany

P +49-(0)30-1208820-89  
F +49-(0)30-1208820-90  
[info@picoquant.com](mailto:info@picoquant.com)  
[www.picoquant.com](http://www.picoquant.com)